

# DRAAGBARE HPGE



# Table of contents

<b>BSI</b> .....	<b>3</b>
MONOLITH Gamma & X-ray HPGe Spectrometer .....	4
HPGe Spectrometer with Lead Shield .....	5
Portable HPGe Gamma- & X-ray Spectrometer .....	6
Robotic Gamma Spectrometer .....	7
Ultra Low-background HPGe Detectors .....	9
<b>PHDS</b> .....	<b>9</b>
GeGI: Imaging HPGe Detector .....	11
NP Radiochemistry Imager .....	12
Fulcrum: HPGe Detector .....	13
Fulcrum-40h: 40% HPGe Detector .....	14
LoPro HPGe Detector .....	15
<b>Overige</b> .....	<b>15</b>
Spectrometer voor snelle gamma-neutronenactivatieanalyse (PGNAA) .....	17



## Partner **BSI**



Baltic Scientific Instruments (BSI) is an OEM manufacturer based in Riga, Latvia, dedicated to the development and production of advanced spectrometric and detection equipment. With decades of experience and roots in the former Research Institute for Radioisotope Apparatus (RNIIRP), BSI provides cutting-edge technologies for nuclear power, environmental monitoring, security, medicine, and scientific research.

The company specializes in HPGe, Si, CdZnTe/CdTe, and scintillation detector systems, known for their accuracy, stability, and performance in demanding analytical environments.

Through continuous innovation, strict quality assurance (ISO 9001:2015), and strong international collaboration, BSI supports customers worldwide in achieving precise and reliable radiation measurement and analysis.

### Product offering


**MONOLITH Gamma & X-ray HPGe Spectrometer**



**HPGe Spectrometer with Lead Shield**



**Portable HPGe Gamma- & X-ray Spectrometer**

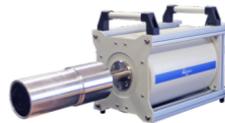

**Robotic Gamma Spectrometer**



**Ultra Low-background HPGe Detectors**





## MONOLITH Gamma & X-ray HPGe Spectrometer

Detection unit Monolith consist from the following integrated components:

- HPGe detector
- Preamplifier
- Autonomous cooling system for the detector based on electrical machinery cooler EMC
- Controller for controlling the operation of EMC
- Fans (2-4) for EMC cooling



### Features

- 10% - 160% efficiency HPGe p-type coaxial detectors are available;
- Energy range from 40 keV to 10 MeV for GCD model;
- Energy range from 3 keV to 10 MeV for GCDX/GCDX-OS models;
- High efficiency of radiation detection;
- High energy rate up to 200000 MeV/sec;
- Excellent peak symmetry;
- Detection of radiation in any spatial orientation depending on cryostat modification;
- Manufacture in a portable cryostat is possible;
- Low background and Ultra - low background materials are available;
- Doesn't require a full thermal cycle after an unexpected shutdown.



## HPGe Spectrometer with Lead Shield

### Application

Coaxial HPGe Detector with Lead Shield is used to measure the specific gamma radiation of radionuclides from various types of environmental objects such as rocks, minerals, sludge, slag, soil, plant, sediment and particulate matter in air and water.

### Features

- Adopting precision gamma-spectrometry methods
- Radionuclide identification and determination of their specific activity
- Low level of instrumental background
- Low threshold for radionuclide detection
- Separate and simultaneous measurement of activity of 100 radionuclides



### DESCRIPTION

Coaxial HPGe Detector with Lead Shield is used to measure the specific gamma radiation of radionuclides from various types of environmental objects such as rocks, minerals, sludge, slag, soil, plant, sediment and particulate matter in air and water.

#### Ordering information

For LN2 cooled dipstick HPGe detection units:

- N100 - 100mm thick lead.
- N150 - 150mm thick lead.

For electrically cooled HPGe detection units (Monolith):

- M100 - 100mm thick lead.
- M150 - 150mm thick lead.

For hybrid cooled dipstick HPGe detection units (Nicole):

- H100 - 100mm thick lead.
- H150 - 150mm thick lead.

← [Back to partner](#)




Stralingsdetectie > Draagbare HPGE

## Portable HPGe Gamma- & X-ray Spectrometer

### Application

Detection, accumulation, processing and analysis of gamma spectra in field and industry conditions where small dimension and weight of spectrometer are important.

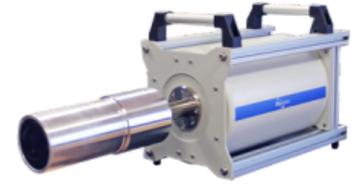

### Features

- Ultra-light cryostat fabrication for minimum gamma absorption;
- Light weight aluminum construction;
- Detection of radiation in any spatial orientation;
- Compact low consuming electronics;
- Available with HPGe coaxial or planar detector;
- Transportation and storage without cooling.

Detection, accumulation, processing and analysis of gamma spectra in field and industry conditions where small dimension and weight of spectrometer are important.



# Robotic Gamma Spectrometer

## Application

The Automated Spectrometer is intended for the detection and analysis of radio nuclides from various types of environmental objects such as rocks, minerals, sludge, slag, soil, plant, sediment and particulate matter in air and water. The spectrometric system is able to determine the composition of a sample based on the photon energy and the activity based on the photon flux. The low-background lead shielding together with the highly pure germanium (HPGe) p-type detector gives precise results even for low activity materials. The fully automated sample changer enables the user to measure more than 40 samples, without having to interact with the Robotic Gamma Spectrometer. This reliable robotic sample changer increases the productivity and reduces the possibility of health risks for the operator.



The Automated Spectrometer is intended for the detection and analysis of radio nuclides from various types of environmental objects such as rocks, minerals, sludge, slag, soil, plant, sediment and particulate matter in air and water. The spectrometric system is able to determine the composition of a sample based on the photon energy and the activity based on the photon flux. The low-background lead shielding together with the highly pure germanium (HPGe) p-type detector gives precise results even for low activity materials.

The fully automated sample changer enables the user to measure more than 40 samples, without having to interact with the Robotic Gamma Spectrometer. This reliable robotic sample changer increases the productivity and reduces the possibility of health risks for the operator.

## Automatic Sample Changer

The seven-axis robot handles a payload of up to 3kg and with, practically, unlimited reach, the robot is able to carry out a series of operations using flexible rather than hard automated solutions. In addition to a horizontal reach, the robot has the ability to reach below its base. Furthermore, the robot has a very compact turning radius, which is enabled by the robots symmetric architecture, without offset on axis 2. This ensures the robot can be mounted close to other equipment.

Basic characteristics of Robot arm:

- Seven-axis manipulator
- Machine vision
- Payload: 3 kg
- Reach: per request
- Fastest 7-axis robot
- Accuracy:  $\pm 0.01$  mm
- Weight: 25 kg
- IP30 protected

- All motors and cablings enclosed
- Compact controller
- Sample holder tool for vessels with diameter in range 40 - 110 mm.
- Barcode Reader and Writer

To assure the correct processing of all data during the measurement and analysis process, the samples are marked by using a barcode printer that is connected to the workstation. Here all necessary information about the sample is stored in a database. Using the bar code reader, the information stored in the database is retrieved for each sample before the measurement process is started. This fully automated process delivers all necessary information for the measurement and analysis process.



# Ultra Low-background HPGe Detectors

## Application

Ultra low-background HPGe detectors are widely used in underground laboratories for determination of radionuclides activities in environmental or industrial samples at  $\mu\text{Bq/kg}$  levels and in scientific experiments such as investigation of magnetic moment of neutrino, dark matter search, etc.



## Design

- Task related design (U-type, vertical, down-looking or portable cryostat)
- Remote not cooled part of preamplifier
- Zeolite is placed near not cooled part of preamplifier in order to be outside measuring chamber

## Cryostat materials

- Certified materials with low radiation impurities
- Ultrapure aluminium-silicon alloy with U + Th content < 1, 0.5 or 0.2 ppb for detector holder and endcap
- Freshly produced electrolytic copper for coldfinger and pedestal
- Tested on radiopurity selected stainless steel screws and sapphire insulators

## Technology

- Transportation of HPGe crystal and cryostat materials by surface freight
- Minimization of fabrication time (location of materials above ground)
- Assembly in a cleanroom
- Cleaning and passivation of copper surfaces
- Storage of crystal and cryostat materials in a container made from materials effectively slowing down and absorbing neutrons (water and Cd)

## Design features

- Fabrication of large volume HPGe detectors without bulletization
- Front end electronics made on low-background Teflon substrate
- Passive screen between front end electronics and HPGe crystal made from Pb with Bi-210 radioactivity < 0.1 Bq/kg
- Double-crystal HPGe detector design
- Multi-crystal HPGe detector design



## Partner **PHDS**



PHDS Co. is gespecialiseerd in de ontwikkeling van draagbare High-Purity Germanium (HPGe) gammastralingsdetectoren. Deze detectoren bieden hoge-resolutie spectroscopie en beeldvormingsmogelijkheden voor toepassingen in nucleaire beveiliging, noodhulp en wetenschappelijk onderzoek. Hun instrumenten zijn ontworpen voor nauwkeurige isotopenidentificatie en -kwantificering in formaten die direct inzetbaar zijn.

### Product offering

**GeGI: Imaging HPGe Detector**



**NP Radiochemistry Imager**



**Fulcrum: HPGe Detector**



**Fulcrum-40h: 40% HPGe Detector**



**LoPro HPGe Detector**



← Back to partner



Stralingsdetectie > Draagbare HPGE

## GeGI: Imaging HPGe Detector

- HPGe gamma-ray spectroscopy and isotope identification
- Real-time exposure rate calculation
- Fully capable of imaging Special Nuclear Materials ( $^{235}\text{U}$  and  $^{239}\text{Pu}$ )
- Fast cool-down to operating temperature
- Long-life mechanical cooler
- Compact and hand-portable
- Hot-swappable batteries for continuous field operation
- User-friendly tablet operation
- Reachback File: ANSI N42.42 format





## NP Radiochemistry Imager

The NP Imager is specifically designed to measure the dynamics of radiochemical separation processes in real time. Developed under a Small Business Innovation Research (SBIR) grant from the Department of Energy Office of Nuclear Physics (NP), the NP Imager focuses on the unique needs of radiochemists and technicians separating radioisotopes for radio-pharmaceuticals and other applications.



The NP Imager monitors the locations and distributions of multiple isotopes over time, allowing the radiochemist and technician to monitor separation processes in real time. In the process shown here, NP Imager measured Lu-177 and Yb-175 separation on a column over the course of 3 ½ hours using 10-minute exposures. Note that it's clear when the Yb-175 has been removed from the column, allowing the radiochemist to know when to collect the desired Lu-177.

### Applications

- Radiopharmaceutical Process Monitoring
- Radiochemical Separation Research and Development
- Nuclear Physics Isotope Production
- Nuclear Materials Management
- Waste Management
- Decontamination and Decommissioning



## Fulcrum: HPGe Detector



### Overview:

The Fulcrum is a next-generation HPGe gamma ray detector engineered for rapid, accurate isotope identification in a compact, field-ready design. Weighing just 8 lbs, it is the lightest and most portable detector of its kind, making it ideal for mobile and emergency response scenarios. Equipped with the user-friendly PHDS OMNI software, the Fulcrum offers real-time exposure rate calculation and intuitive, color-coded isotope identification—all accessible through a simple touchscreen interface. With fast cool-down, wireless capability, and extended field operation features, the Fulcrum sets a new standard for high-performance gamma spectroscopy in any environment.

### Features:

- HPGe gamma-ray spectroscopy and isotope identification
- Ultra-fast cool-down to operating temperature (2 hours)
- Most compact and lightweight HPGe detector (8 lbs)
- Relative efficiency: 12% at 1333 keV
- Real-time exposure rate calculation
- Compact, hand-portable design
- User-friendly Android app interface
- Wireless capability for remote operation
- Long-life mechanical cooler
- Bridge battery option for uninterrupted field use
- User-defined timed data acquisition with auto file save
- Reachback file output in ANSI N42.42 format



← [Back to partner](#)



Stralingsdetectie > Draagbare HPGE

## Fulcrum-40h: 40% HPGe Detector

- HPGe gamma-ray spectroscopy and isotope identification
- Relative efficiency of 40% (at 1332 keV)
- Fast cool-down to operating temperature (5 hours)
- Long-life mechanical cooler
- ISOTAC activity calculator
- Configurable with 1 or 2 onboard batteries and optional neutron detector
- Real-time exposure rate calculation
- User-specified timed data acquisition and auto file save
- Reachback File: ANSI N42.42 format





# LoPro HPGe Detector

The LoPro is a lightweight, low-profile HPGe gamma ray detector used by U.S. military operators around the world. The LoPro provides unmatched field reliability and gamma-ray spectroscopy for mission-critical applications. The unique form factor allows inconspicuous measurements for missions in which discretion is a must, while not sacrificing the spectroscopic performance expected from an HPGe detector.

## Features

HPGe gamma-ray spectroscopy and isotope identification	Fast cool-down to operating temperature (3 hours)	Compact and hand-portable (8 lbs)
Integrated GADRAS isotope identification	Relative efficiency: 12% (at 1333 keV)	Real-time exposure rate calculation
User-friendly Android app, including TAK compatibility	Reachback File: ANSI N42.42 format	Long-life mechanical cooler
Long battery life (up to 15 hours per battery)	Wireless capable for remote operation	User-specified timed data acquisition and auto file save





## Partner **Overige**

---

### Product offering

**Spectrometer voor  
snelle gamma-  
neutronenactivatiean-  
alyse (PGNAA)**





# Spectrometer voor snelle gamma-neutronenactivatieanalyse (PGNAA)

## Toepassing

- De PGNAA-spectrometer is bedoeld voor:
- Analyse van de radionuclidesamenstelling van monsters die sporen van radionucliden bevatten;
- Zoeken naar vervalproducten van radionucliden van antropogene of natuurlijke oorsprong;
- Hoge registratie-efficiëntie en lage detectielimieten worden bereikt met;
- Vermindering van natuurlijke achtergrondniveaus door het gebruik van materialen die zijn gezuiverd van radionucliden;
- Verhoging van het piekcontrast door actieve onderdrukking van Compton-verstrooiingsquanta.

De PGNAA-spectrometer is toepasbaar in:

- Mijnbouw en minerale exploratie: voor nauwkeurige en snelle elementaire analyse van erts en mineralen.
- Milieumonitoring: detecteren en kwantificeren van verontreinigende stoffen in bodem en water.
- Landbouw: analyseren van de bodemsamenstelling om het gebruik van meststoffen en de gewasopbrengsten te optimaliseren.
- Petrochemische industrie: monitoren van de elementaire samenstelling van ruwe olie en derivaten daarvan.
- Geologisch onderzoek: het bestuderen van gesteente- en sedimentmonsters voor wetenschappelijk onderzoek.
- Farmaceutische industrie: het waarborgen van de zuiverheid en samenstelling van grondstoffen en eindproducten.
- Academische en onderzoeksinstituten: het uitvoeren van geavanceerd onderzoek op verschillende wetenschappelijke gebieden.

